# Characterization of miRNA expression in hESC lines using NCode™ SYBR GreenER qRTPCR


**Mark Landers**  
**Epigenetics**  
**Invitrogen**

## Presentation Overview

- **Brief overview of the biology**
- **NCode™ SYBR GreenER miRNA qRTPCR**
  - Platform
  - qRTPCR kit and performance
- **miRNA profiling of hESC cell lines**
  - miRNA expression profiles in hESC
  - Relative Quantitation of miRNA levels in hESC lines

# NC-RNA: The Key to Eukaryotic Complexity?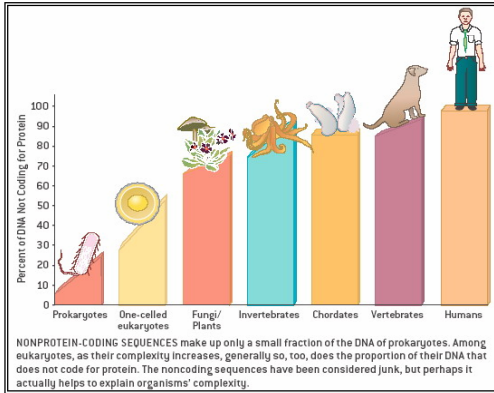

Organism	Percent of Transcriptional Output	
	Protein Coding RNA	Non Coding RNA
<i>E. coli</i>	84	16
<i>S. cerevisiae</i>	71	29
<i>C. elegans</i>	27	73
<i>D. melanogaster</i>	13	87
<i>H. sapiens</i>	2	98


## Possible roles in:

- Cancer
- Neurological disorders
- Host-pathogen interactions

## Data suggesting role in diverse mechanisms:


- RNAi
- DNA methylation
- Chromatin structure/remodeling

# Coding vs. Non-coding RNA

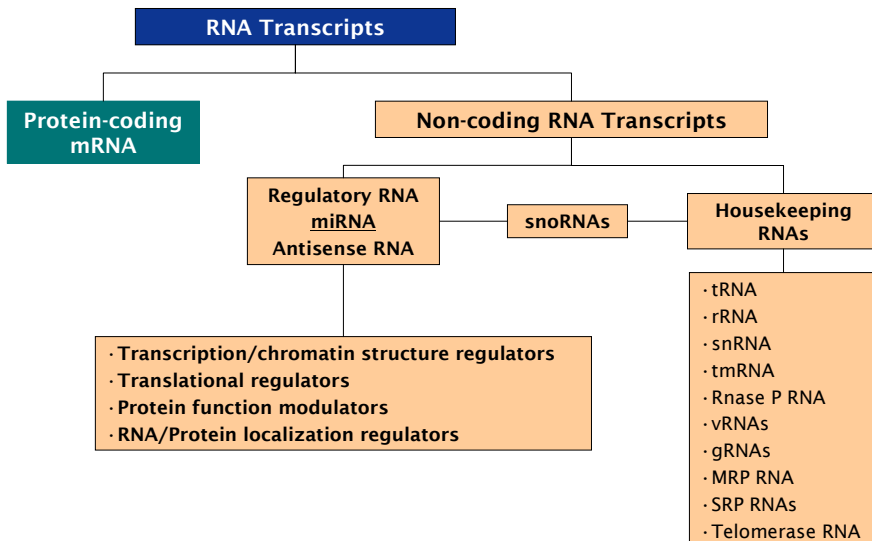NC-RNAs compose majority of transcription in complex genomes

# miRNA Biogenesis and Action
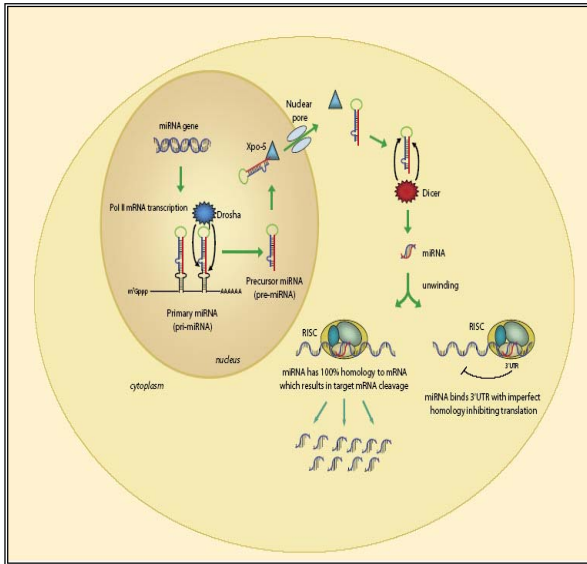
## I. Evolutionarily conserved

- > ~500-1000+ in humans
- > Recent evidence suggests far more could exist...

## II. 20-24 nt in length

## III. Interaction with mRNA target suppressing protein expression

- > Perfect-mRNA degradation
- > Imperfect-translational suppression


## IV. Influence most human protein-coding genes

- > 1/3 mRNAs targeted by miRNAs
- > Remainder of mRNAs "on the run"

5


# miRNA Profiling Workflow

## miRNA profiling workflow overview

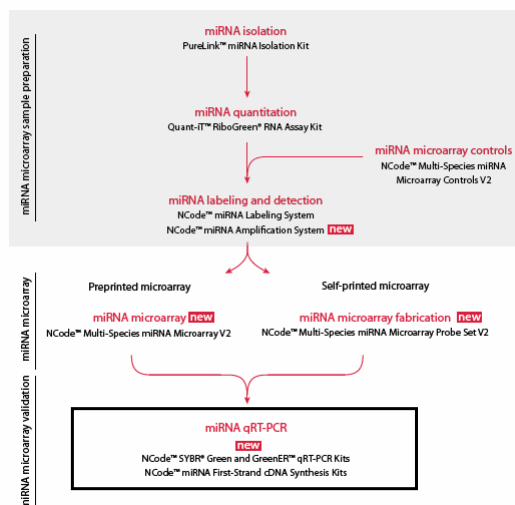6


## NCode™ SYBR® Green miRNA qRT-PCR Kits
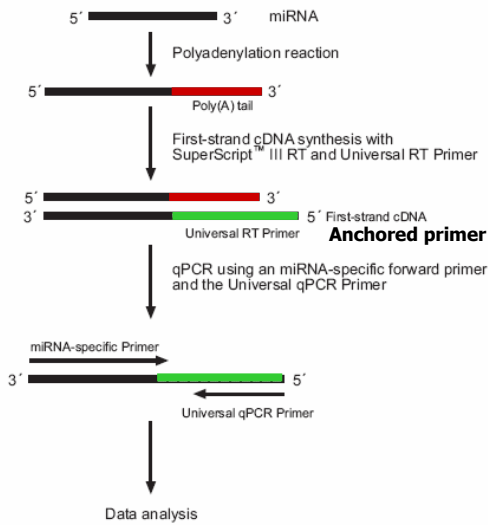
### 1. Flexible system

- Only 1 primer needed
- No special probes
- Quick assay design
- Permits analysis of other small RNAs


### 2. Inexpensive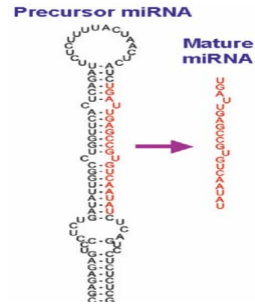

7

invitrogen  
Invitrogen Proprietary & Confidential

## NCode™ SYBR® Green miRNA qRT-PCR Kit

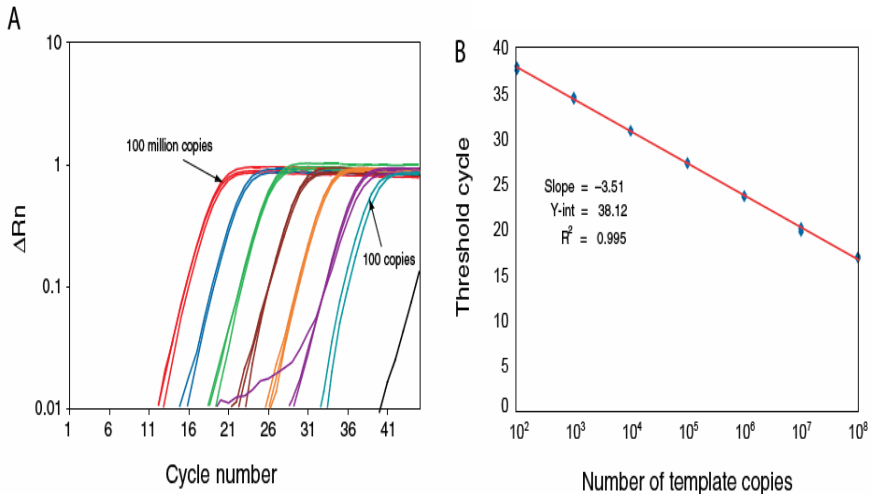
Accurate quantitation over a broad dynamic range of input sample

8

invitrogen  
Invitrogen Proprietary & Confidential

## NCode™ SYBR® Green miRNA qRT-PCR Kit

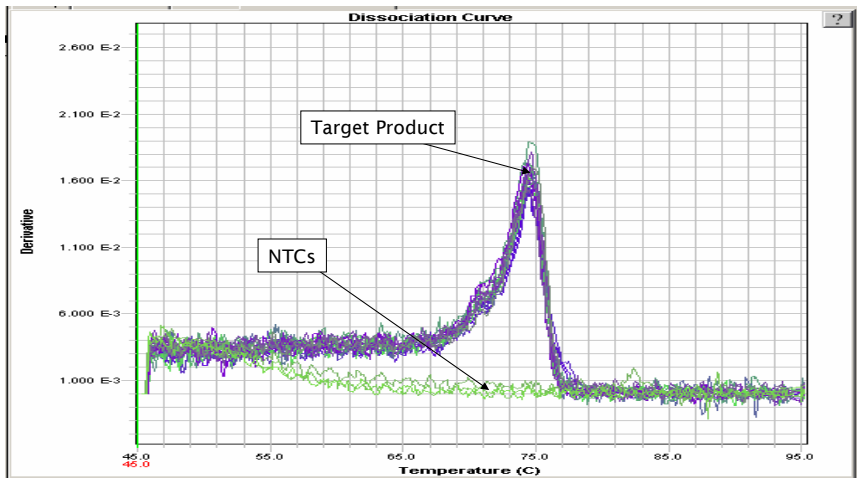
Specific amplification of miRNA targets

9

invitrogen  
Invitrogen Proprietary & Confidential

## Quantification of snoRNAs with NCode™ qRT-PCR

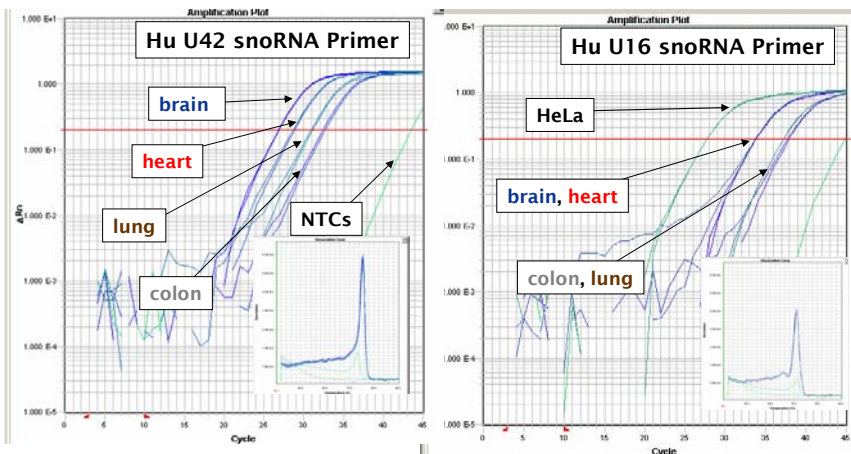
Quantify a range of small RNAs using the NCode™ qPCR system

10

invitrogen  
Invitrogen Proprietary & Confidential

## Housekeeping Genes as Reference Controls

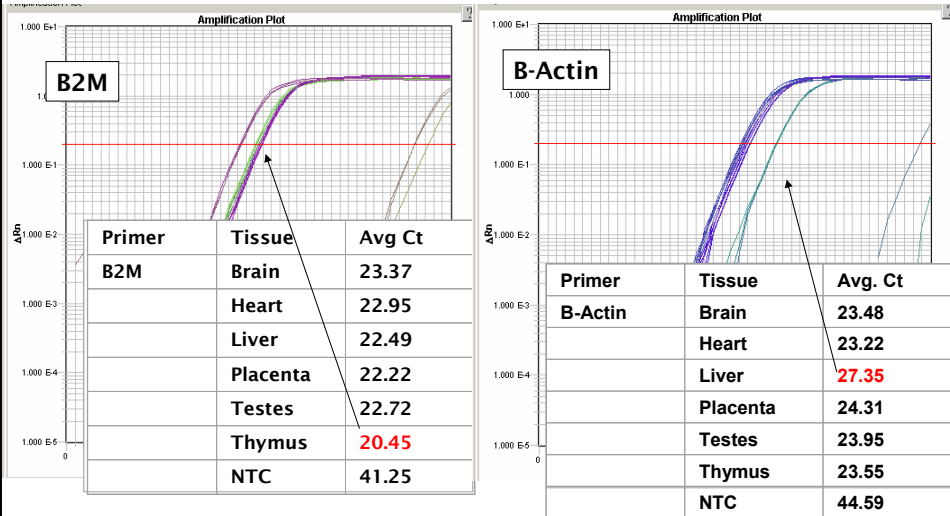
Examine housekeeping gene variation across all potential sample types

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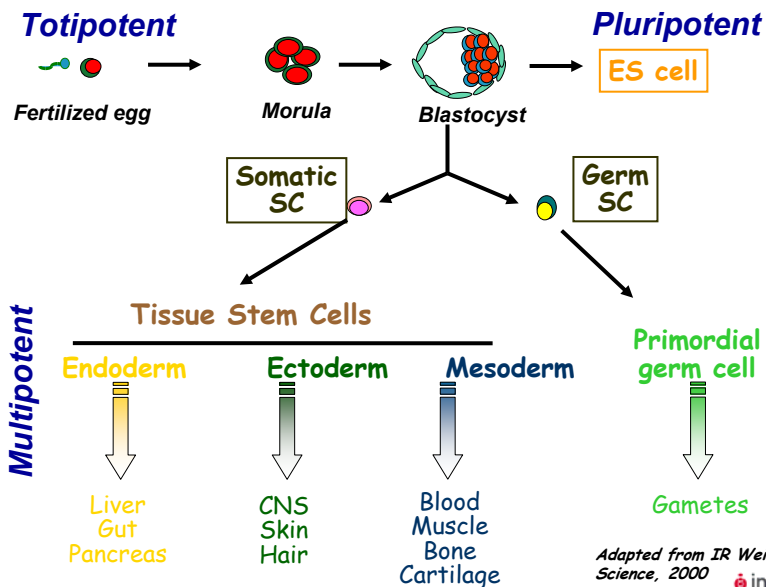

## qRT-PCR: miRNA expression in hESC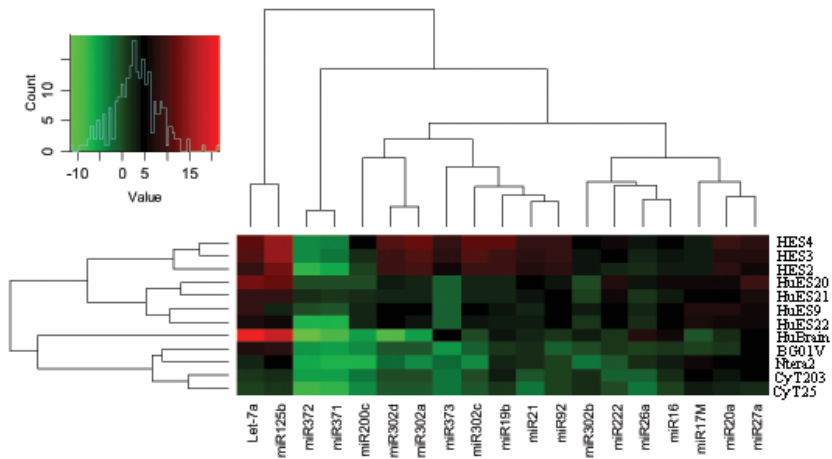

14


## qRTPCR: miRNA expression in ESC vs Differentiated cells

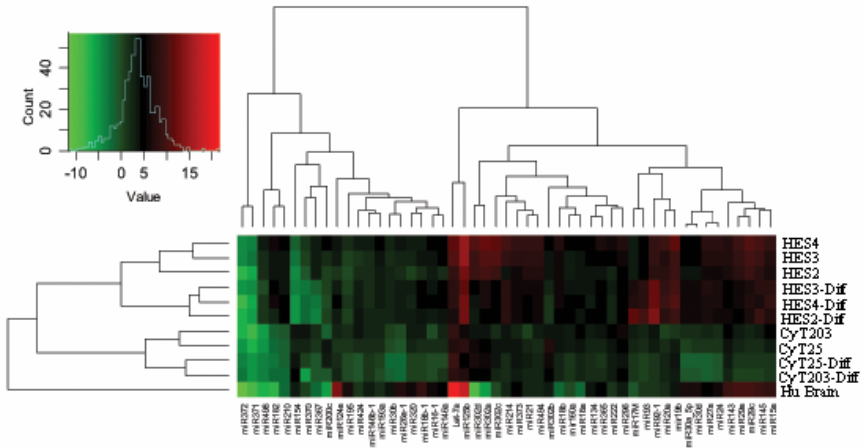15

invitrogen  
Invitrogen Proprietary & Confidential

## Relative miR copy number: hESC vs. EB & Adult brain

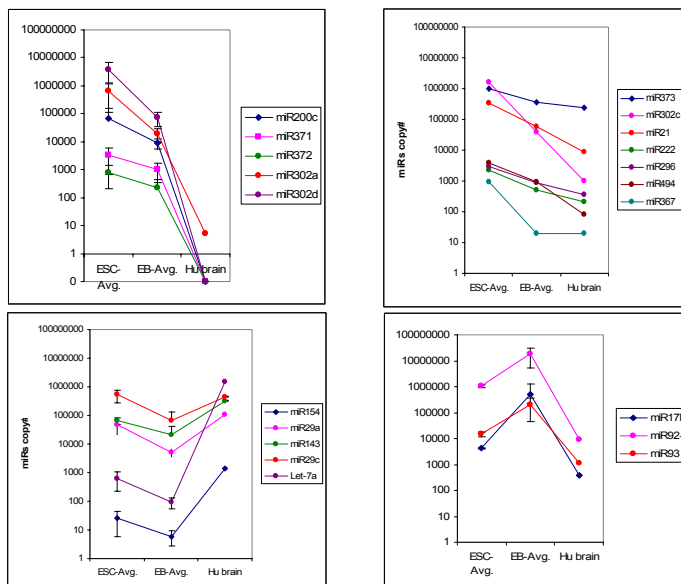16

invitrogen  
Invitrogen Proprietary & Confidential

## miRNAs specific to hESC or EB based on Ncode qRT-PCR profiles of five independent hESC cell lines

miRNA	ESC	EB	Adult Tissue
<i>miR200c, 371, 372, 302a, 320d</i>	<b>High</b>	Low	<i>Absent</i>
<i>miR373, 302c, 21, 222, 296, 494, 367</i>	<b>High</b>	Low	<i>Lower</i>
<i>miR154, 29a, 143, 29c, Let7a</i>	<b>High</b>	Low	<b>High</b>
<i>miR17M, 92, 93</i>	Low	<b>Higher</b>	Low
<i>miR16, 134, 145, 30a_5p</i>	Same	Same	Same

## ACKNOWLEDGEMENTS

### Invitrogen R&D

#### *Stem Cells and Regenerative Med.*

Uma Lamshimpathy  
 Brad Love  
 Bhaskar Thyagarajan  
 Ying Liu  
 Soojung Shin  
 Haipeng Xu  
 Shayne Boucher  
 Kate Wagner  
 Jonathan D. Chesnut  
 Mahendra S. Rao

#### *Epigenetics Group*

Chris Adams  
 Lisa Wenrich

### Collaborators

#### *Rutgers University, NJ*

Ronald P. Hart  
 Loyal Goff  
 Rebecca Jorsten  
  
 Melissa Carpenter, Novocell  
 Ralph Graichen, ES Cell int.

## NCode™ qPCR Kits

- NCode™ SYBR® GreenER™ miRNA qRT-PCR Kit
  - Advanced SYBR® GreenER™ technology
  - 10 poly(A) rxns, 20 cDNA rxns, 100 qPCR rxns
- NCode™ SYBR® Green miRNA qRT-PCR Kit
  - 10 poly(A) rxns, 20 cDNA rxns, 100 qPCR rxns
- NCode™ miRNA First-Strand cDNA Synthesis Kit
  - 10 reaction kit
  - 50 reaction kit